# **SAFETY, STORAGE AND SETUP INFORMATION**

## **PRODUCT SUPPORT SERVICE AND REPAIR**

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## SAFETY, STORAGE AND SETUP INFORMATION

- 1) **General Information:** the following general precautions for handling must be observed during all phases of operation and service of CAEN equipment. Failure to comply with these precautions or with specific warnings elsewhere in the product manual could result in personal injury or damage to the equipment. The precautions listed below represent warnings of certain dangers of which CAEN is aware. The user of the product, should follow these warnings and all other safety precautions necessary for the safe operation of the equipment in the operating environment.

**CAEN shall not be liable for user's failure to comply with these requirements.**

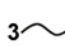
- 2) **Conventions Used in This paragraph:**

**CAUTION indicates a hazard that could damage the equipment. Such damage may invalidate the warranty.**

- 3) **Safety Terms and Symbols on the Product:** these terms may appear on the product:

- **DANGER:** indicates a hazardous situation which, if not avoided, could result in serious injury or death.
- **WARNING:** indicates a hazardous situation which, if not avoided, could result in death or serious injury.
- **CAUTION:** indicates a situation or condition that, if not avoided, could cause physical injury or damage the product and/or its environment.

The following symbols may appear on the product:

		
Caution, refers to the product manual	Caution, risk of electric shock	Productive Conductor Terminal
		
Earth (Ground) Terminal	Alternating Current	Three-Phase Alternating Current

- 4) **Unpacking and Inspection:**

**CAUTION:** observe precautions for handling electrostatic sensitive devices.

- Inspect containers for damage during shipment. Report any damage to the freight carrier for possible insurance claims.
- Check that all the components received match those listed on the enclosed packing list. (CAEN cannot accept responsibility for missing items unless we are notified promptly of any discrepancies.)
- Open shipping containers; be careful not to damage any contents.
- Inspect contents and report any damage. The inspection should confirm that there is no exterior damage to the unit such as broken knobs or connectors and that the front panel and display face are not scratched or cracked. Keep all packing material until the inspection has been completed.
- If damage is detected, file a claim with carrier immediately and notify CAEN service.
- If equipment is to be installed later, place equipment in original shipping container and store in a safe place until ready to install.

- If equipment must be returned for any reason, carefully repack equipment in the original shipping container with original packing materials if possible. Please contact CAEN service.
- **CAUTION:** some equipment **MUST** be shipped with the original packing case (this is indicated in the manual and on the packaging). **Failure to comply may cause damage to the equipment and may void the warranty.** In this case, please save and store all packaging materials.

**5) ESD precautions:** CAEN equipment may be classified as ESD sensitive item<sup>1</sup>.

**CAUTION:** during long term or temporary storage the items shall be covered by ESD protective bags. Handling of the items without ESD protective covering shall be performed only into approved ESD Protected Area (EPAs); any operator shall be properly grounded (e.g., by wrist-straps) before handling the item. Avoid handling operations if the relative humidity is below 30% or take extra ESD safety precautions (such as redundant ground clips) if such handling is unavoidable.

<sup>1</sup> Excluded all the devices provided with independent case or chassis. Devices that consist of a circuit board with no case or chassis (e.g. VME, CAMAC, NIM modules or HV/LV Power supply boards) comply with the EMC directive if installed in a CE marked system.

**6) Storage Condition<sup>2</sup>:** CAEN equipment shall be handled and stored in a clean, dust free environment under the following conditions:

	Minimum	Maximum
<b>Temperature</b>	- 30° C	+ 80° C
<b>RH (non-condensing)</b>	5 %	90 %

<sup>2</sup>Typical ranges, for more details see the product manual.

**7) Handling:** CAEN equipment shall be handled only with lint-free gloves (e.g. nitrile); never touch the item with bare, ungloved hands.

**CAUTION:** do not subject the item to undue shock or vibrations; do not bump, drop or slide shipping containers. Do not leave items or shipping containers unsupervised in areas where untrained personnel may mishandle the items.

**8) Instruction for Cleaning:** if required, the equipment may be cleaned with isopropyl alcohol or deionised water and air dried. Clean the exterior of the product only. Do not apply cleaner directly to the items or allow liquids to enter or spill on the product.

**Cleaning the Touchscreen (if present)**

In order to clean the touchscreen, wipe the screen with a towelette designed for cleaning monitors or with a clean cloth moistened with water. Do not use sprays or aerosols directly on the screen; the liquid may seep into the housing and damage a component. Never use solvents or flammable liquids on the screen.

**Cleaning the air vents (if present)**

It is recommended to clean the air vents (if present) occasionally on all vented sides of the board. Lint, dust, and other foreign matter can block the vents and limit the airflow. Be sure to unplug the board before cleaning the air vents and follow the general cleaning safety precautions.

**CAUTION:** CAEN recommends cleaning the device using the following precautions:

- Never use solvents or flammable solutions to clean the board.
- Never immerse any parts in water or cleaning solutions; apply any liquids to a clean cloth and then use the cloth on the component.
- Always unplug the board when cleaning with liquids or damp cloths.
- Always unplug the board before cleaning the air vents.
- Wear safety glasses equipped with side shields when cleaning the board

**9) Transport:**

**CAUTION:** CAEN equipment can be safely transported in their original shipping packages.

**CAUTION:** some equipment **MUST** be transported with the original packing case (this is indicated in the manual and on the packaging). **Failure to comply may cause damage to the equipment and may void the warranty.**

**10) Installation and Operation:** only qualified personnel should perform installation, operation, and maintenance procedure.

**CAUTION:** Periodically verify the proper functioning of the equipment before use to avoid misinterpretations or incorrect readings that could compromise safety.

**CAUTION:** Installation must ensure proper connection to the power supply, protective grounding, the use of any required external protective devices, adequate ventilation, and compliance with noise emission limits.

**CAUTION:** For equipment with a mass greater than 18 kg, use the integrated lifting handles.

**CAUTION:** Maintenance must be performed only by qualified personnel; after each intervention, carry out the appropriate safety checks.

**11) Safety Precautions:** keep extreme caution when an electrical shock hazard is present. Lethal voltages may be present on cable connector jacks or test fixtures. Warning label indicates HV sources which are hazardous to personnel; during test or integration all exposed HV connections shall be insulated or protected. Before operating an instrument, make sure the line cord is connected to a properly grounded power receptacle. Inspect the connecting cables, test leads, and jumpers for possible wear, cracks, or breaks before each use. When installing equipment, where access to the main power cord is restricted, such as rack mounting, a separate main input power disconnect device must be provided, in close proximity to the equipment and within easy reach of the operator. It is recommended to disconnect ac line from the entire test system and discharge any capacitors before:

- Connecting or disconnecting cables or jumpers.
- Installing or removing switching cards.
- Removing cover and making internal changes, such as installing or removing jumpers, if required.
- **CAUTION:** The instrument and accessories must be used in accordance with its specifications and operating instructions, or the safety of the equipment may be impaired.
- **CAUTION:** Use only the supplied accessories.
- **CAUTION:** Do not exceed the maximum signal levels of the instruments and accessories, as defined in the specifications and operating information.
- **CAUTION:** When fuses are used in a product, replace with same type and rating for continued protection against fire hazard.

**12) Ventilation Requirements**

**CAUTION:** all power supplied CAEN NIM, VME and CAMAC modules must be used only with crates with forced cooling air flow.

CAEN modules without power supply, such as Attenuators and Delay units, do not require crates with forced cooling air flow.

**CAUTION:** rack bound CAEN HV Power Supply Systems must be installed with a minimum 10 cm space both above and below the system.

**13) Medical Use:** CAEN equipment is not designed with components and testing procedures that would ensure a level of reliability suitable for use in treatment and diagnosis of humans. These items are not intended to be a substitute for any form of established process or equipment used to monitor or safeguard human health and safety in medical treatment.

#### 14) Operating Environmental Requirements<sup>3</sup>

	Minimum	Maximum
<b>Temperature</b>	0° C	+ 40° C
<b>RH (non-condensing)</b>	10 %	90 %
<b>Environmental</b>	Indoor use	
<b>Altitude</b>	<2000m	
<b>Pollution degree</b>	2	
<b>Overvoltage category</b>	II	
<b>EMC environment</b>	Commercial and light industrial	
<b>IP degree</b>	IPX0, not for wet location	

<sup>3</sup> Typical ranges, for more details see the product manual.

#### 15) Installing Devices with No Case or Chassis: (e.g., VME, CAMAC, NIM modules or HV/LV Power supply boards):

**CAUTION:** CAEN strongly recommends to use an antistatic wrist strap and a conductive foam pad when installing these items. After removing the devices from the system or its protective antistatic wrapper, place the component flat on a grounded, static-free surface (and in the case of a board, component side up). Do not slide the component over any surface. Do not stack units on top of each other.

**CAUTION:** crate power should be turned off during insertion or removal of modules<sup>4</sup> to avoid possible damage caused by **momentary** misalignment of contacts.

**CAUTION:** the safety of any system that incorporates the device falls under the responsibility of the assembler of the system

<sup>4</sup> Excluded modules supporting live insertion/extraction into/from the crate.

**CAUTION:** to insert a module in the relative crate follow these steps:

- Choose a convenient I/O station.
- Hold the board firmly with both hands (if HV/LV Power supply boards).
- Align the PCB back side with one crate free slot rails.
- Start sliding gently the board along the rails.
- Pay attention to closer boards: insert the module carefully to avoid contacting boards to the right or left.
- Ensure that correct alignment is kept.
- Continue sliding the board holding it firmly.
- Ensure that sliding is smooth.
- Plug the board into Crate backplane connectors with gentle finger pressure.
- Ensure the correct backplane connectors are mating.
- Complete installation by screw-locking the board front panel on the crate.

#### 16) Repairs: may be executed by CAEN service only.

#### 17) EMI considerations: CAEN equipment has been designed in compliance with the harmonized standards concerning Electromagnetic compatibility. Devices meet requirements concerning EMC. Specific test was performed in normal use condition in order determine the effective EM effects. Adhering to the following rules will help EMI.

- Power supply cable routing. The power supply cable is usually susceptible to conduct EMI inside the device and coupling into power supply cord. It is recommended to keep the power supply cord far from SIP/SOP cables and fixed in place with cable ties, if necessary.

- SIP/SOP connector shield (if available). Sometimes the I/O connector realize poor contact between metal housing and connector plate. It is recommended to check electrical continuity.
- Minimize the differential mode loop of signal wires by keeping the wires close together. This will minimize the magnetic field and reduce radiated EMI.
- Minimize the common mode capacitance of wires by keeping the length as short as possible.
- If specified, connect the chassis earth connection properly.
- If applicable, the devices are designed to run with spread spectrum clocking enabled: it is recommended because often produces low EM emissions.
- If not necessary, disable unused clocks. Clock signals that have no load can have high level of ringing that can lead to EMI problems. For the same reason it is recommended to disable unused HV channels or input acquisition channels.

**CAUTION:** If the customers modify the hardware/software/firmware factory configuration, the EMC compliance will not be guaranteed.

**18) Efficient Energy Use – General Guidelines:** to promote efficient use of electricity and reduce environmental impact, it is recommended to:

- Turn off the device when not in use.
- Optimize battery usage by avoiding partial or unnecessary charging (if applicable).
- Charge the device only when needed based on battery level, using appropriate chargers (if applicable).
- Disconnect the device from the power source once charging is complete or usage has ended.
- Keep the device in good working conditions to ensure consistent energy efficiency.

**19) Device Decommissioning:** After its intended service, it is recommended to perform the following actions:

- Detach all the signal/input/output cable.
- Wrap the device in its protective packaging.
- Insert the device in its packaging (if present).

**CAUTION:** the device shall be stored only at the declared environmental condition, otherwise performance and safety will not be guaranteed.

## 20) VME Crates / VME Digitizers compatibility

**CAUTION:** do not use the CAEN V/VX1740, V/VX1751, V/VX1730, V/VX1743, V/VX2740, V/VX2745, and VX2730 digitizers with VME8001, VME8002, VME8004, and VME8004A crates: overheat may damage the modules.

The following table shows the compatibility of all CAEN VME CRATES with all digitizers.

		Digitizers																							
		VX2730	VX2751	V2740	VX2740	V2745	VX2745	V1720	VX1720	V1724	VX1724	V1725S	VX1725S	V1730S	VX1730S	V1740	VX1740S	V1742	VX1742	V1743	VX1743	V1751	VX1751	V1761	VX1761
VME64X Crates	VME8004X	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
	VME8008X	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
	VME8100-VME64X <sup>5</sup>	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
	VME8200	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
	μ-crate	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓	✓
VME64 Crates	NV8020A			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8004B			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8008B			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8010			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8011			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8100-VME <sup>6</sup>			✓		✓		✓		✓		✓		✓		✓		✓		✓		✓		✓	
	VME8001							✓		✓		✓						✓						✓	
	VME8002 <sup>7</sup>							✓		✓		✓						✓						✓	
	VME8004 <sup>7</sup>							✓		✓		✓						✓						✓	
	VME8004A <sup>7</sup>							✓		✓		✓						✓						✓	

<sup>5</sup> Models of the VME8100 series equipped with VME64X backplane.

<sup>6</sup> Models of the VME8100 series equipped with VME64 backplane.

<sup>7</sup> Obsolete models.

For the latest version of this table, refer to the following link:

<https://www.caen.it/faq/caen-vme-crates-digitizers-compatibility-table/>

## 21) Regulatory Compliance

- EMC: CE 2014/30/EU Electromagnetic Compatibility Directive.
- Safety: CE 2014/35/EU Low Voltage Directive.
- RED: CE 2014/53/EU Radio Equipment Directive (if applicable).

## 22) Use of Rechargeable Batteries: If the device is equipped with rechargeable batteries, follow the instructions below:

- Use only the battery supplied with the device.
- Recharge according to the manufacturer's instructions..
- Recharge using the dedicated charger and in a well-ventilated area. Do not leave the battery charging unattended.
- Do not disassemble, short-circuit, expose to heat, fire, or excessive moisture.
- Replace the battery only if specified and strictly follow safety instructions.
- Do not use or recharge the battery if damaged, swollen, or overheated.
- Immediately stop using it if the device shows abnormal behavior.
- Dispose of the battery at the end of its service life only through authorized collection centers, in accordance with applicable regulations.

## 23) Preamplifiers: Some preamplifiers operate with detectors that require high-voltage bias. For these units, the following precautions apply:

**CAUTION:** During normal operation, a potentially hazardous high-voltage bias is applied to the detector through the preamplifier. Installation, operation, and maintenance of this unit must be carried out only by qualified personnel.

The preamplifier bias circuit has a very long time constant, meaning that the circuitry can remain charged at high voltage for an extended period. If adequate caution is not exercised, this residual voltage may cause personal injury due to electrical shock. Please observe the following precautions:

- Completely discharge the detector bias circuit by switching off the bias supply before connecting any cable to the input/detector connector.
- When using a variable power supply, reduce the output voltage to zero and wait at least 30–60 seconds. The bias circuitry will then discharge itself through the bias supply output.

**CAUTION:** Do not connect the input/detector connector to exposed circuitry. Connect the preamplifier to a Detector/Power Supply properly grounded to safety earth.

**24) High Voltage Power Supplies:** This family of devices generates high voltages; therefore, specific precautions are required.

**CAUTION:** do not power a load outside its specified range.

**CAUTION:** do not connect or disconnect HV cables while energized or connected to a voltage source.

**CAUTION:** do not operate this product with covers or panels removed.

**CAUTION:** do not operate this product in wet or damp conditions.

**CAUTION:** do not operate this product in an explosive atmosphere.

**CAUTION:** do not operate with suspected failures. if you suspect that this product is damaged, contact technical support.

**CAUTION:** the HV output generates extremely hazardous voltages with potentially lethal currents. Do not connect or disconnect HV cables when the power switch is ON. Always switch off the power and discharge any residual voltage, before connecting or disconnecting HV cables.

**25) Disposal:** The disposal of the equipment must be managed in accordance with Directive 2012/19 / EU on waste electrical and electronic equipment (WEEE).

The crossed bin symbol indicates that the device shall not be disposed with regular residual waste.

## PRODUCT SUPPORT SERVICE AND REPAIR

**For technical assistance on hardware and software, please follow the following procedure**


- 1) **MyCAEN Account:** A MyCAEN account is required for ask support about CAEN products and register a firmware/software license. Making an account is totally free and it is possible to upgrade to MyCAEN+ anytime.

### **New user registration**

- Go to registration page and fill in the registration form with your data.
- Validate your email address by clicking on the confirm link and set a password.
- Log in to CAEN website to activate your account.

### **Do I have a MyCAEN+ account?**

If you are already registered and need to know if you have a MyCAEN+ accounts, you will be able to check it by logging in to CAEN website, go to your dashboard and see the title of the section that could be MyCAEN or MyCAEN+.

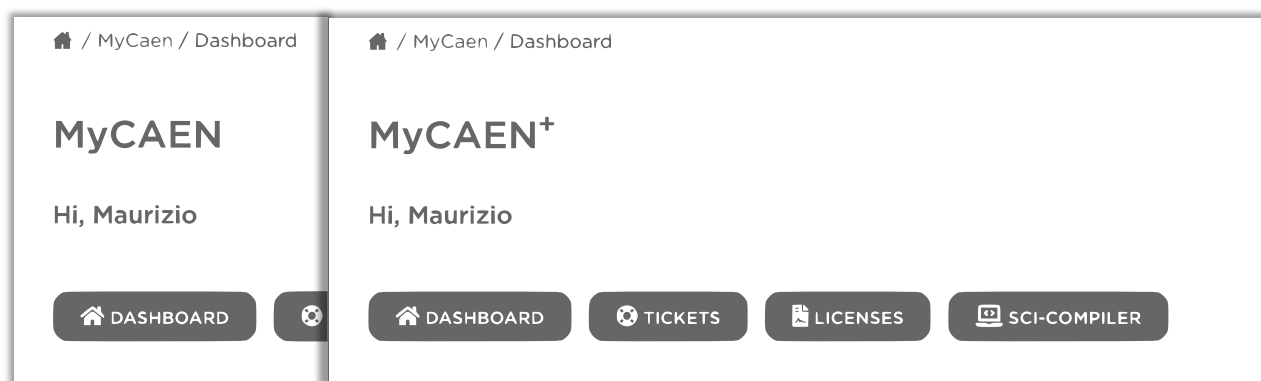### **Upgrade to MyCAEN+**

If you are already registered on CAEN website and you have a MyCAEN account, you will just need to update to MyCAEN+ by following the instructions below:

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.


- Go to "Edit Profile" and on "Profile Information" update your Institutional email, country/region and Company/institute fields.
- Click on "Save Profile".

Known institutes are automatically certified by CAEN website. New institutes or generic email addresses need to be validated by CAEN staff. This operation can take up to 48 hours.

2) **Register a Device.** The product registration is required for access to support services.

## Add a device

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button on the top right.
- On the menu area click on “Dashboard”, then “Add Device” button.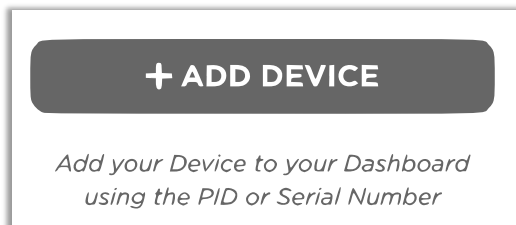


- Select an identification method, PID or S/N. Please refer to chapter “Find your S/N or PID” to discover how to find your board identification number.

The form is titled 'Select an identification method' with a magnifying glass icon. It has two tabs: 'PID' (selected) and 'S/N'. Below the tabs are input fields for 'PID\*' (containing '12345') and 'Model Name'. A 'SEARCH' button is to the right. Below the fields is a note: 'Insert your PID or S/N and click the Search button for device correspondences.' There is an 'Ordering Option' field. At the bottom, there is a checkbox for 'I have read the information about the Privacy Policy and I agree the use of my personal data.' and an 'Add new device' button.

- Click on “Search” button and then “This is my device.”

The result is titled 'Products corresponding to this SN/PID'. It shows a device icon, the model 'NIM8302', and the description 'NIM 5U Compact crate, 10 slot, 150W'. Below this, it shows 'Ordering Option: WNIM8302150W' and 'SN/PID: 12345'. A 'This is my device' button is on the right.

- The fields will be automatically filled by the system. If all information are correct and the Privacy Policy is accepted, it will be possible to click on “Add new device”.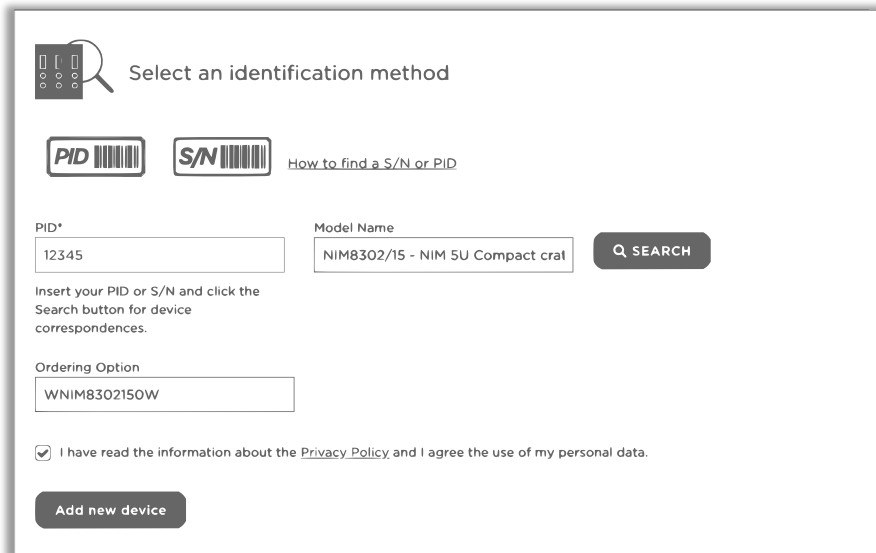

Select an identification method

**PID** **S/N** [How to find a S/N or PID](#)

PID\*  Model Name

Insert your PID or S/N and click the Search button for device correspondences.


Ordering Option

☒ I have read the information about the [Privacy Policy](#) and I agree the use of my personal data.

If the device cannot be found, please double check your identification number; if the search gives no results instead, it will still be possible to create a support ticket without registering the device: the model name and the Serial Number or PID can be specified in the message.

### Find your Serial Number (S/N) or PID

The product information label contains your Serial Number (S/N) or Product Identifier (PID). A Serial Number is typically composed of two or more numbers, while a PID consists of at least five characters. This label is located either on the back of your CAEN module, on the front, or on the underside of the right side of the board.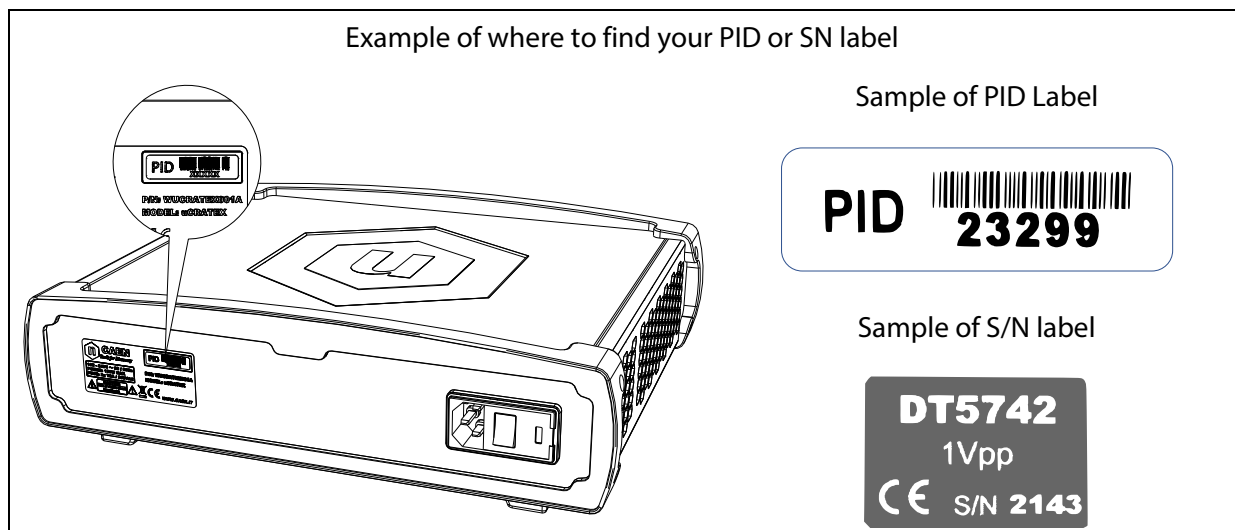


For a detailed guide on how to locate the Product Identifier (PID), please refer to the following link: <https://www.caen.it/find-your-pid>

### Download device list (.csv)

It is possible to download the device list in .csv format, by the following:


- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Dashboard", then "Download devices list" written on the right.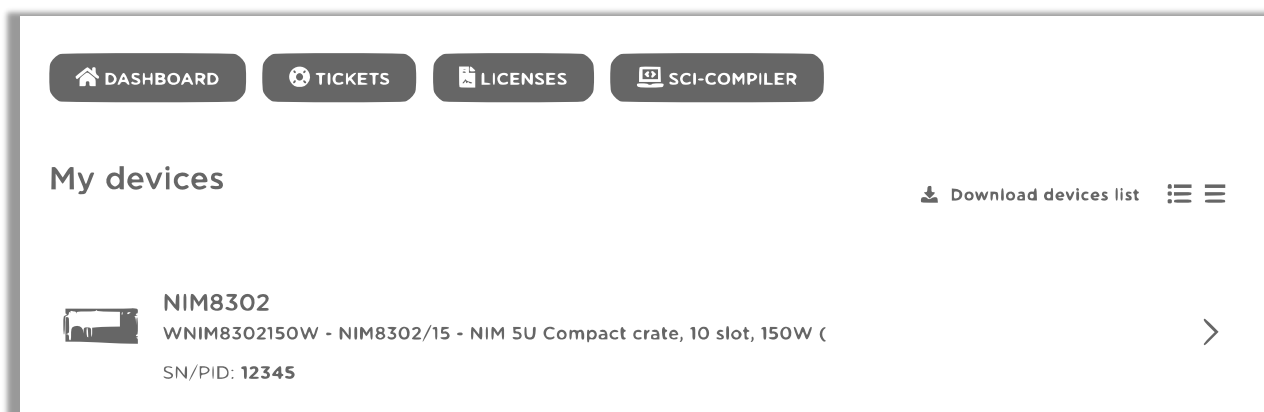

### 3) Support Ticketing

#### Create a Ticket

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Tickets", then "Create Ticket" button.

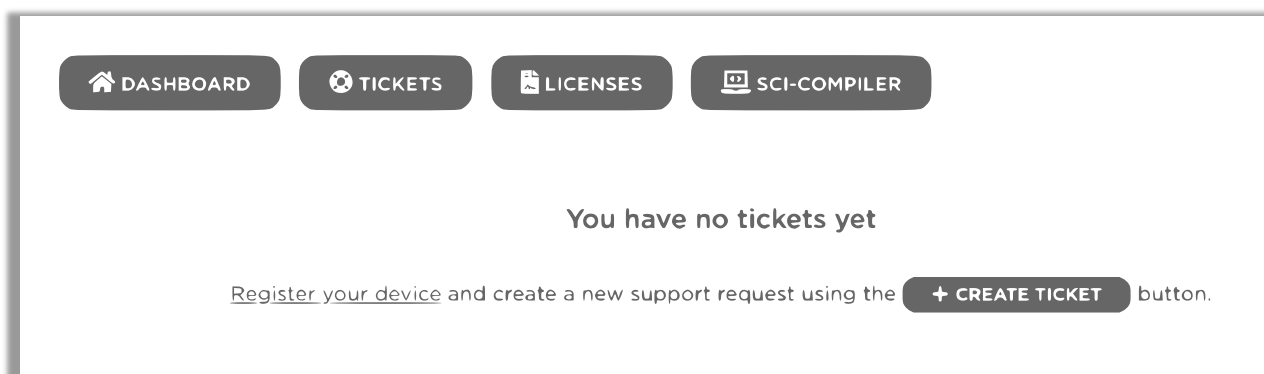- Fill all the required field being careful to select your hardware device.

When the CAEN team reply to your tickets you will receive a notification email with a direct link to the relative ticket.

#### Reply to a Ticket

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "Tickets", then click on the tickets you want to reply.
- Fill the form at the bottom of the page.

When the CAEN team replies to your tickets you will receive a notification email with a direct link to the relative ticket.

#### RMA – Return Merchandise authorization

- Create a Ticket describing your hardware failure.
- If your problem is evaluated by CAEN team as not solvable by the user, the RMA form will be activated, and a notification email will be sent to the user.
- Click on the link received by email and then click on "Return Device" to fill the form with all return information.

Depending on where you are located, the product must be shipped directly to CAEN S.p.A. headquarters or to your Country Dealer. To find the proper address please visit the page at

<https://www.caen.it/sales-network/>

4) **License.** The user can register the software license.

**Web License**

- Log in to CAEN website and go to your dashboard by clicking on MyCAEN button at the right top.
- On the menu area click on "License", then "Register License".
- Complete the wizard procedure by selecting the module in your possession.

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